## **VALENCIACOLLEGE**

## Chemistry

## Lab Technique 23: Fan-Folding a Flat Filter Paper (Video starting at 1:04.)

Fold flat filter paper in half three times to create quadrants of equal size.	Triater raper (video starting at 1.04.)
Open once.	
Then make two more folds by bringing the outer edges into the centerfold.	
Open it completely.	
To create fan-folds, take three fold lines. Push the center fold down and bring the two outer folds together. Repeat all the way around the filter paper until it is completely pleated.	Up Down
Open it keeping pleats.	